

Long term storage of *Alternaria*

June 2014

Evenhuis, Wageningen

Trudy van den Bosch and Marieke Förch, Plant Research International.

Long term storage of *Alternaria* species

Put 1 ml sterile 15% DMSO (diluted in tap water) in a sterile 2 ml cryovial. Add a few pieces of the *Alternaria* infected agar plate (1-2 week old plate) or *Alternaria* spores. Put the cryovial in a Styrofoam box overnight at -80°C. Put after 24 hours the cryovial in a liquid nitrogen canister.